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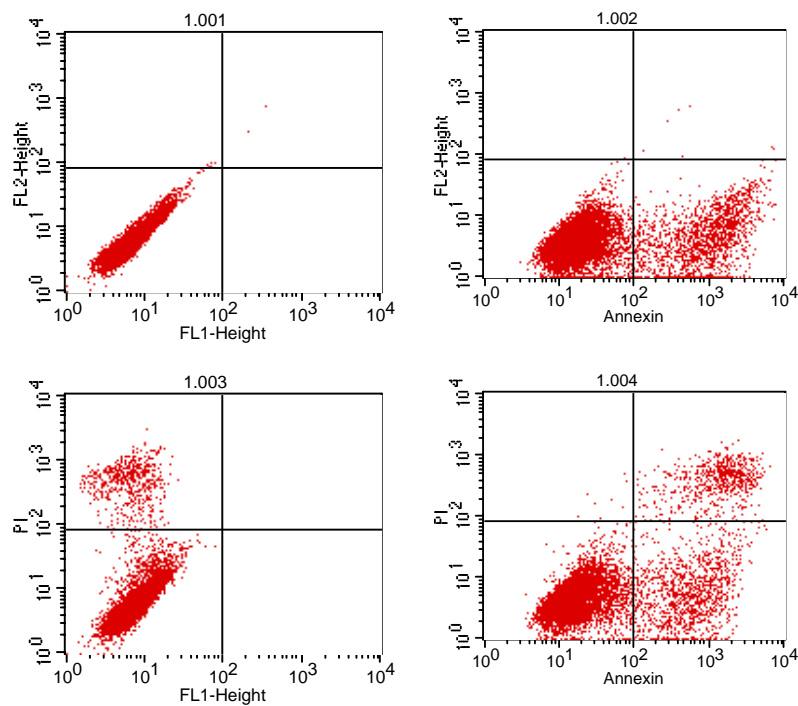
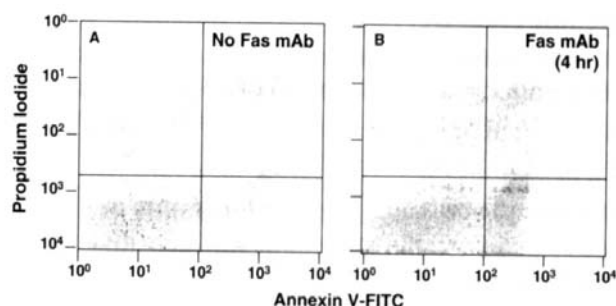
Figure 12. Flow Cytometric Analysis of Apoptotic Cells Using Annexin V-FITC. Jurkat T cells were left untreated or were treated for 4 hr with anti-human Fas mAb, clone DX2 (Cat. No. 33450D) and Protein G. Cells were incubated with Annexin V-FITC in a buffer containing PI and analyzed by flow cytometry. (A) Untreated cells were primarily Annexin V-FITC and PI negative (lower left quadrant, panel A), indicating that the cells were viable and not undergoing apoptosis. (B) After treatment with DX2 and Protein G, a significant number of cells are Annexin V-FITC positive and PI negative (lower right quadrant, panel B), indicating that the cells are in an early stage of apoptosis and are still viable. In addition, a population of cells has progressed to a later stage of apoptosis and are stained with both PI and Annexin V, indicating that the cells are no longer viable.

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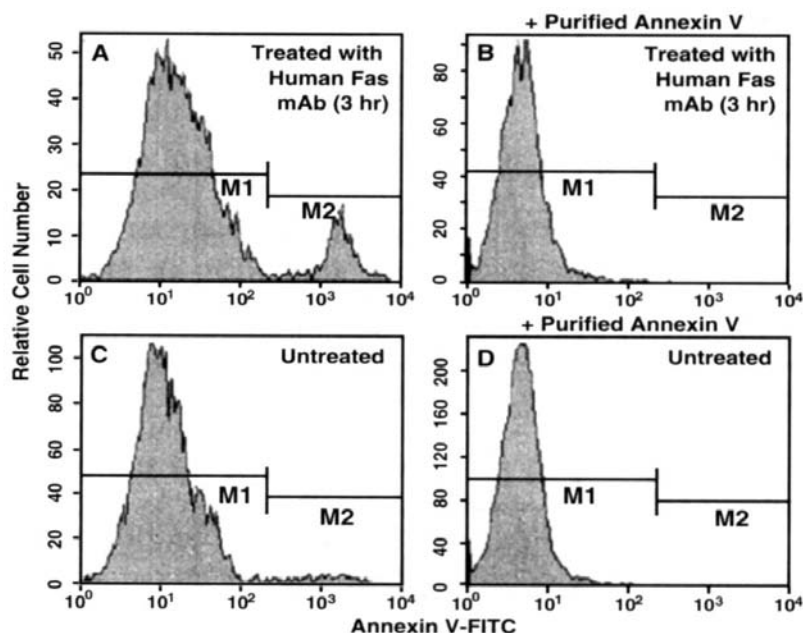
Figure 11. Flow Cytometric Analysis of Annexin V Staining and Blocking. Jurkat T cells were induced to undergo apoptosis by treatment with anti-Fas antibody, clone DX2 (Cat. No. 33450D) and Protein G for 3 hr (A, B). Other cultures were left untreated (C,D). Cells were incubated with Annexin V-FITC alone (A,C) or with recombinant Annexin V to block Annexin V binding sites prior to adding Annexin V-FITC (B, D). After a 3 hr treatment with DX2, a population of cells was Annexin V-FITC positive (A, M2 gate). Annexin V-FITC staining was blocked when cells were first incubated with purified recombinant Annexin V (B). As expected, the cell population that was not treated with DX2 were primarily Annexin V negative (C). The small number of Annexin V positive cells in the untreated population likely represents a basal level of apoptosis. This was blocked when cells were first incubated with purified recombinant Annexin V (D). The M1 and M2 gates demarcate Annexin V-FITC negative and positive populations, respectively.

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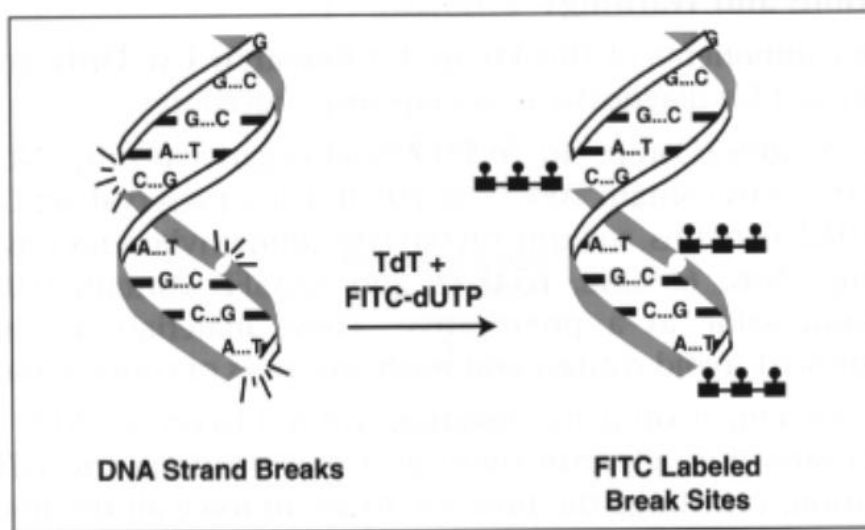
Schematic Representation of APO-DIRECT™ Labeling. The enzyme terminal deoxynucleotidyl-transferase (TdT) catalyzes a template independent addition of FITC-labeled deoxyuridine triphosphates (FITC-dUTP) to the 3'-hydroxyl ends of double- and single-stranded DNA. DNA break sites are identified by the FITC-labeled dUTP's.

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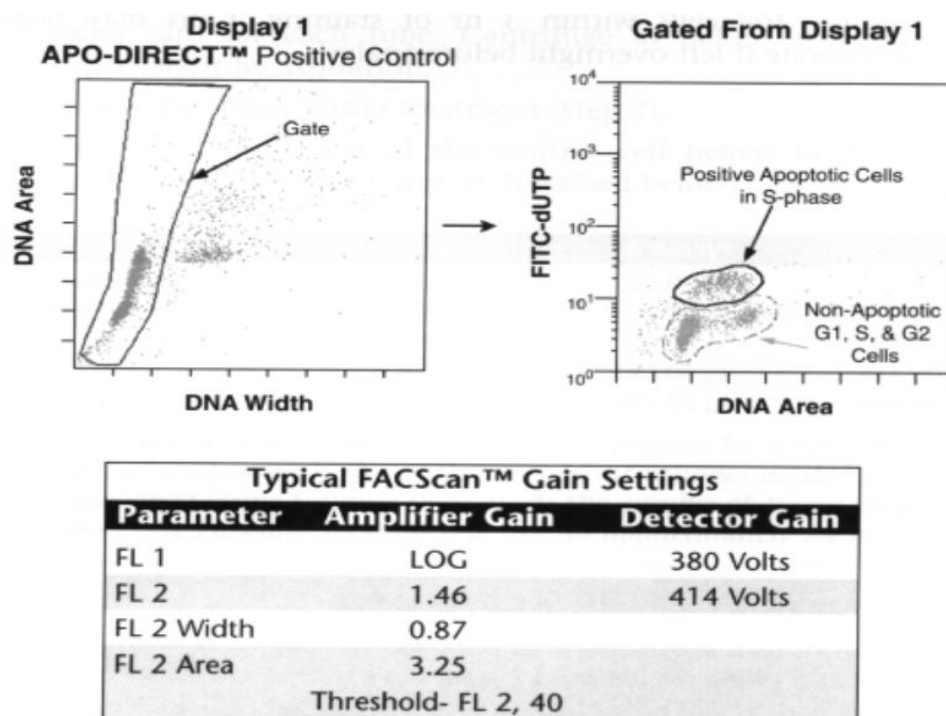
Figure 17. Flow Cytometer Setup for Becton Dickinson Hardware. Cells are labeled with both PI (DNA) and FITC (Apoptotic Cells). Display 1: Non-clumped cells are gated. Gated from Display 1: A circle is drawn around cells that stain positive with FITC. An example of FACScan™ gain settings is shown. Note that in this cell population, apoptotic cells are in S-phase. However, apoptosis may also occur in other phases of the cell cycle.

5. 阴性质控细胞 Negative Control Cells 和阳性质控细胞 Positive Control Cells:

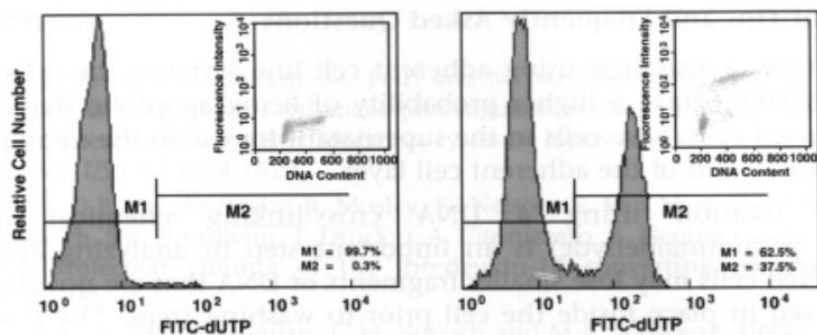
Figure 18. Flow Cytometry Data of APO-DIRECT™ Negative and Positive Control Cells. Typical results using the positive and negative control cells are shown in this figure. Negative control (left panel) or positive control cells (right panel) are treated with FITC-dUTP in the presence of TdT enzyme which incorporates FITC-dUTP into the 3'-OH-DNA ends found in apoptotic cells. Cells are then stained with propidium iodide and analyzed on a Becton Dickinson FACScan™. The presence of apoptotic cells in the right hand panel is demonstrated by increased fluorescence intensity (M2 gate). Propidium iodide staining (panel insets) provides additional cell cycle information about each population, indicating that the majority of apoptotic cells are in S-G2 phases of the cell cycle.

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APC BrdU Flow Kit

Cat. No. 552598 (50 tests) Cat. No. 557892 (4×50 tests)

Component	Storage condition	No. of vials
Fluorochrome -anti-BrdU Antibody	4°C	1
BD Cytofix /Cytoperm Buffer	4°C	1
BD Perm/Wash Buffer (10×)	4°C	2
BD Cytoperm Plus Buffer	4°C	1
BrdU (10 mg/ml)	-80°C	5
DNase	-80°C	5
7-AAD	4°C	1

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每管50 μl 染色50 样本(10^6 cells/样本). FITC BrdU Flow Kit(Cat. No. 559619) 包含50 μl FITC anti-BrdU抗体。APC BrdU Flow Kit (Cat. No.

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Sample Data using the BD Pharmingen FITC BrdU Flow Kit Protocol

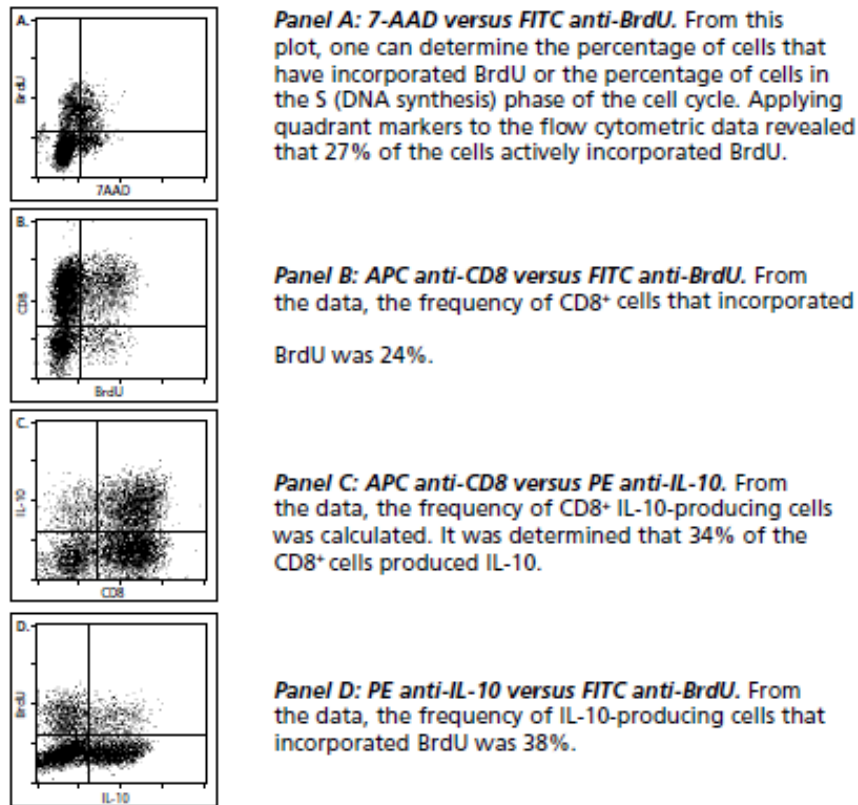
Figure 1. Multicolor flow cytometric analysis of stimulated mouse cells that synthesized DNA and/or produced IL-10. Spleen cells from a BALB/c mouse were primed *in vitro* and restimulated with PMA and ionomycin in the presence of a protein transport inhibitor (to promote intracellular cytokine accumulation). During the final 45 minutes of culture, the cells were labeled with 10 μ M BrdU. The cells were then harvested and stained with FITC anti-BrdU (FL1), PE anti-IL-10 (FL2), 7-AAD (FL3), and APC anti-CD8 (FL4). The panels depict two-color staining patterns generated from the reanalyzed flow cytometric data for these cells. A great deal of information can be obtained for a given cell population stained using this multiparameter staining technique. (see panel descriptions)

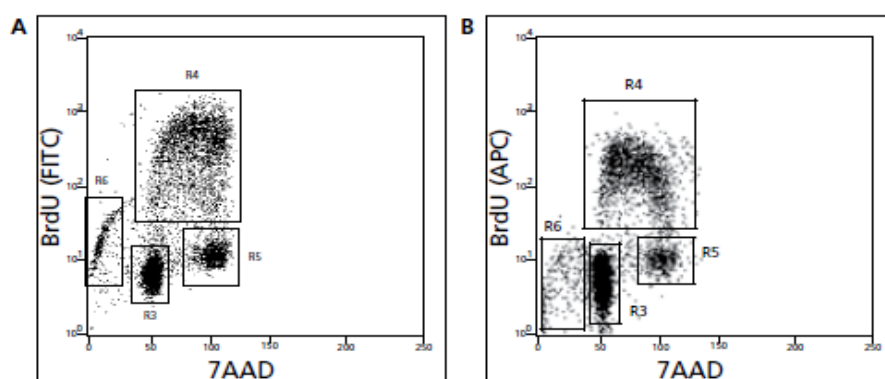
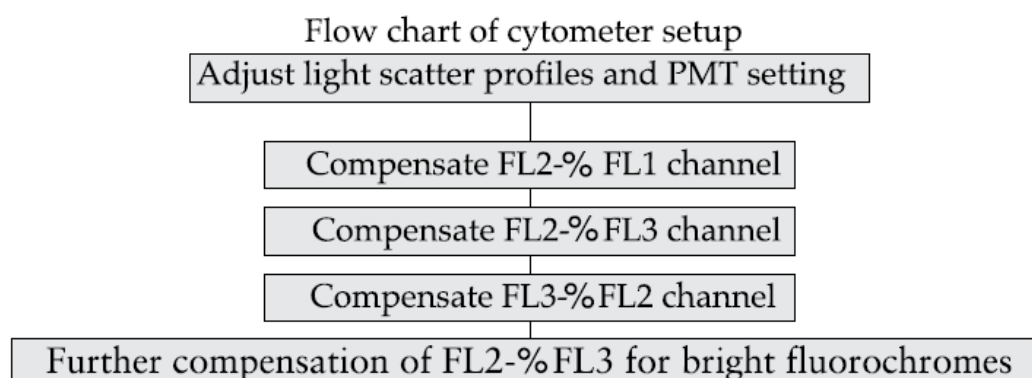
Figure 3. Region gates for the quantitative cell cycle analysis of populations that have been stained for incorporated BrdU and total DNA levels. **Panel A:** The measurement of cell incorporated BrdU (with FITC anti-BrdU) and total DNA content (with 7-AAD) in D10.G4.1 cells. The D10.G4.1 cells were cultured with 10 μ M BrdU for 30 minutes. The cell cycle positions and active DNA synthetic activities of cells can be determined by analyzing the correlated expression of total DNA and incorporated BrdU levels. As shown by the region gates applied to the 7-AAD versus BrdU dot plot, flow cytometric analysis of cells stained with the reagents provided in the BrdU Flow Kit allowed for the discrimination of cell subsets that were apoptotic (defined as sub-G0/G1, R6, 5.6% of cells) or resided in G0/G1 (R3, 38.6%), S (R4, 38.6%), or G2 + M (R5, 14.4%) phases of the cell cycle and had recently synthesized DNA. **Panel B:** The measurement of cells-incorporated BrdU (with APC anti-BrdU) and total DNA content (with 7-AAD). Human PBMCs were stimulated with immobilized anti-human CD3 antibody, clone HIT3a, 10ug/ml for plate coating, (Cat. No. 555336), soluble anti-human CD28 antibody, clone CD28.2 at 2ug/ml, (Cat. No. 555725), recombinant human IL-2 at 10ng/ml (Cat. No. 554603) and recombinant human IL-4 at 20ng/ml (Cat. No. 554605) for 2 days. The cells were then washed and subsequently expanded for 3 days in culture with medium containing recombinant IL-2 and IL-4. Finally the cells were harvested and restimulated for 4 hours with PMA (Sigma, Cat. No. P-8139 5ng/ml) and ionomycin (Sigma, Cat. No. I-0634 500ng/ml). 20uM BrdU was added for the final hour. Region gates are as in figure 3A with region 6 being apoptotic (3.31%), region 4 is S phase(23.5%), region 3 is G0/G1(64.3%), and region 5 is G2+M (6.1%).

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